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0.5ml holder

Manufacturer	Sumitomo Bakelite	As one
Product Name	Proteosave SS	PCR tube
Type	MS-4201X MS-4205M MS-4255M	2-8086-02
Material	Sumiron proteosave	Polypropylene
Compatibility	OK	OK

1.5ml holder

Manufacturer	Sumitomo Bakelite	As one
Product Name	Proteosave SS	Low adsorption tube
Type	MS-4202X MS-4215M MS-4265M	2-3971-02
Material	Sumiron proteosave	Polypropylene
Compatibility	OK	OK

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15ml holder

Manufacturer	Falcon	Sumitomo Bakelite	As one
Product Name	Falcon tube (PP)	Proteosave SS	Centrifuging tube
Type	352096 352097 352196	MS-52150	2-8089-11(ECK-15ML-R)
Material	Polypropylene	Sumiron proteosave	Polypropylene
Compatibility	OK	OK	OK

Manufacturer	NUNC	Kartell
Product Name	Centrifuging tube	Conical tube
Type	366036 339650	84000 84001 84006 84007 84008 84010
Material	Polypropylene	Polypropylene
Compatibility	OK	OK

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30ml holder for Careful Cleaning(NotForSample)

Manufacturer	SARSTEDT
Product Name	Centrifuge tube with Cap 30ml
Type	60.544 60.544S(Sterilized) 60.545(Sterilized)
Material	Polypropylene
Compatibility	OK

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5ml holder

Manufacturer	Falcon	BM Equipment
Product Name	Falcon tube (PS)	Test tubes 12x75mm 4.0ml
Type	352003 352008 352054 352058 352235	BMT-1275-PP
Material	Polystyrene	Polypropylene
Compatibility	OK	OK

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Manufacturer	NUNC
Product Name	Centrifuging tube
Type	366036
Material	Polypropylene
Compatibility	OK

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8 well strip plate holder

Manufacturer	Thermo Scientific	Sumitomo Bakelite	Corning International
Product Name	Piece 8-well strip plates	Separate plate 8 FM	PCR tube strips 8 well
Type	15031	MS-8508M	9102
Material	Polystyrene	Polystyrene	Polystyrene
Compatibility	OK	OK	OK

96 well plate holder

Manufacturer	Falcon	Falcon	Falcon
Product Name	6 well micro plate	12 well micro plate	24 well micro plate
Type	351146	351143	351147
Material	Polystyrene	Polystyrene	Polystyrene
Compatibility	OK	OK	OK

Manufacturer	Falcon	Falcon
Product Name	48 well micro plate	96 well micro plate
Type	351178	351172
Material	Polystyrene	Polystyrene
Compatibility	OK	OK

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Manufacturer	Corning
Product Name	384 Well Clear Flat Bottom
Type	#3702
Material	Polystyrene
Compatibility	OK

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96 PCR plate holder

Manufacturer	Roche Diagnostics	Roche Diagnostics
Product Name	LightCycler 8-tube strips	LightCycler multiwell plate 96
Type	6327672	4729692
Material	Polystyrene	Polystyrene
Compatibility	OK	OK ※製造停止 (230925確認)

→メーカー動作確認済み (230925追記)
LightCycler® 8-Tube Strips (white)
(Roche Diagnostics 06612601001)

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Manufacturer	Roche Diagnostics	Roche Diagnostics	Bio-Rad
Product Name	LightCycler multiwell plate 384, White	LightCycler multiwell plate 384, Clear	Hard-Shell 384-Well 480 PCR Plates
Type	4729749001	05102413001	HSR-4801
Material	Polypropylene	Polystyrene	Polystyrene
Compatibility	OK	OK	OK

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